PHANTOM HIGH-SPEED CAMERA

Miro C321/C321J/ N5/JBox 2.0 USER MANUAL





When it's too fast to see, and too important not to.®







The contents of this manual are subject to change without notification.

PN: ZDOC-64106-MA-0003 Rev 2 Last Updated: August 2022

PHANTOMHIGHSPEE

When it's too fast to see, and too important not to.®

Miro C321/C321J/N5/JBox 2.0 User Manual

1	Introduction	1
2	Connectors	5
3	Quick Start Guide	9
4	Battery	13
5	PCC Phantom Software	15
6	Signals & Programmable I/O	25
7	Miro Junction Box 2.0	29
8	Accessories	33
9	Support	41

CONTENTS

D.COM



VEO 440 & Miro C321J

The Phantom Miro C321, C321J, N5 camera head and Junction Box 2.0 (JBox 2.0) are the building blocks for flexible multiple-camera setups.

- Either Stand-alone or with a system for flexibility: The Miro C321 and the N5 camera head with the Miro N-B Base can be used either individually or as part of a multi-camera system.
- Multi-camera system efficiency: The Miro C321J and N5 with Miro N-JB Base connect to the Junction box with 1 simple, rugged cable. Power, signaling and data all flow through the Junction Box.
- Mix and match with Phantom VEO cameras: The Junction Box 2.0 is compatible with up to 4 VEO cameras as well as up to 6 of all other Miro C cameras and legacy Miros.
- Expandable: The Junction box can be daisy-chained or branched to as many junction boxes as the application needs.

Their small size, rugged body, and flexibility make them perfect for applications including:

- Vehicle safety testing
- Ballistics testing
- Microscopy
- Mechanical analysis

CAMERA CONTROL

Phantom cameras come with PCC software for Ethernet-based setup and control.

IMAGE STORAGE

The C321 and C321J have either 8GB or 16GB of internal memory, and the N5 bases have 8GB of memory. Both cameras and bases have a 240GB internal, non-removable CineFlash to protect data in the event of power loss.

BATTERY BACKUP

The Miro C and N5 bases have an internal, non-removable battery designed to provide back-up power to complete operation and safely save a cine in the event of power loss to the camera. The battery is UNDOT 38.3 qualified.

IMAGE MONITORING & VIDEO OUTPUTS

Phantom Miro C cameras have a 1.5G HD-SDI output on the front of the camera via a mini-DIN connector. The N5 has 1.5G HD-SDI on the base via a standard BNC connector. In each case, a video monitor or view finder can be connected to view the image at the camera. Standard video rates up to 1080 p30 are supported.

ADDITIONAL FEATURES

Set new CSR default: Set a CSR in PCC with customer-selected settings. If settings are not changed, the camera may be powered off and on and the CSR will remain active. Camera must be set at full resolution.

Programmable I/O: Assign and define signal parameters for each Programmable I/O port. Available signals include Strobe, Ready, F-Sync, Pre-Trigger, Event, Memory Gate and several core signals. See Programmable I/O section for more information.

Data Acquisition: External sensor data imported through a compatible National Instruments DAQ model can be embedded in Cine files and visualized in PCC along with the high-speed images.

Image-Based Auto-Trigger (IBAT): Trigger the camera (or a number of connected cameras) from motion detected within the live image.

Exposure Tools: Auto-exposure, Threshold mode for the video output, PCC histogram, and a Zebra pattern overlay in PCC are all tools used to dial in the best exposure without over saturation.

Multi-Cine: Support for up to 63 partitions.

Quiet Fans: When enabled, the fans are stopped during capture to eliminate vibration.

Continuous Recording: Automatically and continuously saves Cines to external storage.

Shutter Off Mode: (also known as Exposure in PIV) turns off the shutter to maximize achievable exposure with minimal frame overhead.

Burst Mode: Generate a precise number of frames with every frame sync pulse.

DHCP: Dynamic Host Configuration Protocol enables the automatic assignment of an IP address when using a DHCP router. This method eliminates the requirement to define the PC's dedicated camera network. Automatic DHCP is an alternative to the secondary user-definable IP network available in all Phantom cameras.

HV Trigger: High Voltage (HV) trigger, when selected in PCC the camera will trigger from a +6-32V pulse.

Detailed information about Phantom cameras, features, and software can be found at: *www.phantomhighspeed.com*



Miro C321J Rear View



Miro C321 Rear View



Miro NJB Top View



Miro NB Top View



Miro C321J Front View



Miro C321 Front View



Fischer connector connects to a Camera Port on a Miro Junction Box (J-Box).





BATTERY RESET

Recessed push-button turns battery power off with use of paper clip or similar object.



Fischer connector connects GigE to a control unit computer / laptop for camera control communication.



Fischer connector connects to a Miro MiniBoB or standard Miro adapter cable.



Fischer power connector connects to standard 40W Miro Power supply.



Capped BNC connector for HD video on N5 bases.

- CXP BNC connector for CXP to N5 camera head.
 - SDI DIN 1.0/2.3 connector for HD video.

C321 AND MIRO N BASE **CONNECTORS**



MIRO JUNCTION BOX

A READY

BNC connector; closed' for ready, 'open' for not ready; covers all cameras connected to this JBox and any JBox up-chain of this one.

B TRIGGER

BNC connector; contact closure causes trigger. BNC connector also accepts TTL pulses (rising / falling edge) for trigger. There is also Trigger on System connector: isolated, ←1V deassert, 3 to 24V assert.



Selectable strobe or FSYNC for synchronization with equipment such as pulsed LED lighting.

D ETHERNET

Ruggedized RJ45 connector connects GigE to acontrol unit computer / laptop for camera control communication.



Same Fischer connector as Camera port. Connects to a Camera Port on second JBox to daisy chain or branch the system for additional cameras.

Fischer connectors; each Miro JBox has six (1-6) Camera ports which carry

be used to daisy chain or branch to another JBox's System port.

Ethernet, power, serial, and Miro camera signals on capture connector. There are two status LEDs for each camera port – Ethernet and camera status. One port can

POWER 20-36VDC Independent power is required for each Junction Box.



QUICK START GUIDE

Vision Research, Inc. | Miro C321/C321J/N5/JBox 2.0 Camera Manual



Phantom cameras are typically controlled with PCC software through a dedicated Ethernet network. PCC is compatible with the 64-bit version of Windows 10 operating systems.

The latest version can be downloaded at: **www.phantomhighspeed.com/PCC**

ASSIGNING A Gb CAMERA NETWORK

Connecting the camera using Gb Ethernet simply requires the computer to detect the IP address range of Phantom cameras.

PCC software includes a **Network Configuration Utility** which will launch automatically after the software is first installed, or it can be later launched from the PCC program itself.

The NetConfig Utility will show all network interfaces currently installed in the PC. Identify the one that will be physically connected to the camera (this process can be done with or without a camera connected at the time).

Click 'Config Adapter', which will then adjust the IP Address and Subnet to the first one available within the 100.100 network that all Phantom cameras are pre-configured to work with. Once connected, the Phantom logo will appear on the lower right with the "Configured for PCC operation" message.

OPTIONAL: MANUALLY ASSIGN Gb CAMERA NETWORK

Alternatively, In the Windows 'Network and Sharing Center,' select the camera network. Change the Ethernet IP settings by selecting 'Properties,' then 'TCP/IPv4' 'Properties,' then select 'Use the following IP address.'

Enter IP address **100.100.100.1** and **Subnet 255.255.0.0**. Other settings should be blank.

POWER ON CAMERA

Connect the camera's dedicated power supply to the Primary DC input or connect via system cable to the Miro JBox. The camera is ready for operation when the red capture light of the Trigger button is illuminated.

LAUNCH PCC SOFTWARE

Double-click the PCC icon located on the desktop. Camera will be recognized immediately if connected and network settings are correct.

SELECT CAMERA FOR USE

In the '**Manager**' tab, select the Phantom camera to be used from the '**Cameras**' group folder. Click the '**Live**' tab.

DEFINE RECORDING PARAMETERS

Click '**Cine Settings**' and define the following parameters by either selecting the value from the pull-down selection list or typing a value into the respective data entry field.

- 1. Set '**Resolution**' to the required Width x Height.
- 2. Choose the required 'Sample Rate,' and 'Exposure Time'
- 3. Set '**Post Trigger**' to zero (0) by moving the '**T**' (Trigger Position) slider to the right, or enter zero (0) into the '**Last**' data entry field.

Cover the lens and click the 'CSR' button to perform a Current Session Reference. A CSR is a black reference for that session. It is required before capturing the first cine and after changing any recording parameters.

With color cameras, perform a White Balance by right-clicking an area of the image that is neutral gray or white, as long as it is not 100 percent saturated.

SET DEFAULT CSR

For customers that rarely change their camera settings, the "**Set new CSR default**" feature provides a convenient alternative to this process. It is available in versions of PCC 3.77 and higher. The camera must operate at full resolution, for example 1920 x 1080 for the C321. The Set Default CSR is set for the frame rate and exposure time being used. That CSR will remain through turning the camera off and on, for as long as that setting is not changed.

Step 1: Enable in PCC:

ook Canera Enable	v Set To: Enable v Hy	physe Group	Look.	Bold -	r.
Show Menus	Enable v 🛛 Expand 1	Neru On Sta	n		
Canera Settings	Controls	Appears	ince	Group L	ock
Cine Settings	Lock Cine	Enable	¥		~
Rash Memory	Set Alt	Enable	v		11
Advanced Settings	Name	Enable	¥		
Salley Control	Description	Enable	~		
Att Force m	Resolution	Enable	¥	Yes v	
Frame Rate Profile	Samuela Rater	Enable	-	Nes w	
Image Based Auto-Trigger	Functione Times	English		Nam or	
Continuous Recording	Expense mile	Englis		Ten	
Camera Ho	Farmer index	English		Max. Lot	
	Expense model	Enable .		198 1	
	LIN COLUMN THE COLUMN	LT-SOLO	*	1088 V	
	C3K options with right click:	race	*	-	
	Low Light	Enable	Y	568 V	w

Step 2: Right click the CSR button in the live tab, and select "Set Default CSR"

Exposure Ind	ex 2000 ~								
EI + Image Processing 2000									
<u>C</u> S Image R	Set new CSR default CSR With Shutter Open CSR With File Save								
-25358	Last: V 0								

FINE-TUNE SETTINGS

After CSR and White Balance are performed, adjust settings, aperture and/or lighting to get a good exposure. A CSR must be performed after any camera settings are adjusted.

'ARM' CAMERA

Click the 'Capture' button to start recording to the camera's internal RAM memory buffer.

TRIGGER

At the end of the action, click the '**Trigger**' button at the bottom of the '**Live**' panel provide an external trigger signal (TTL pulse) via the Trigger connector.

PLAYBACK AND EDIT CINE

Click the '**Play**' tab to Scrub through the timeline or use the '**Video Control**' buttons to locate the first image to be saved.

Click the 'Mark-in' 📧 button.

Locate the last image of the Cine to be saved and then click the 'Mark-Out' 🔟 button.

SAVE TO FLASH OR COMPUTER

Click the large 'Save Cine...' button on the bottom of the 'Play' panel.

- 1. Navigate to the folder where you want to save the Cine file.
- 2. Enter a file name and from the 'Save as Type', select the 'Cine Raw, .cine' file format.
- 3. Click the 'Save' button to begin downloading the Cine file from the camera.

CONFIRM COMPUTER SAVE

Click the '**Open File**' 📂 button. Navigate to the folder and open the saved Cine file. Review the playback by scrubbing through the file and viewing the playback.

BATTERY



OPERATION

The Miro C cameras and N5 Bases have an internal, non-removable battery designed to provide back-up power to complete an operation and safely save a cine in the event of power loss to the camera or base. The following description of battery operation pertains to both Miro C cameras and the N5 bases. Only for simplicity, the description refers only to the camera. The battery is UNDOT 38.3 qualified and has a total capacity of 8.4 volts. When the battery activates, it ready to provide power to the camera in the event DC power is lost, and the LED is Green. The following conditions must be met for battery to activate:

- The camera has DC power and is fully booted.
- The battery is charged to at least 8.0 volts.
- In the Battery Control Section in PCC, the box "Enable Battery" is checked (Default).

The 3 modes of operation are:

Preview Mode (Selectable): This mode provides a 2-minute window to disconnect and reposition cameras during set-up without needing to re-boot the cameras. The battery activates when the camera is placed in Preview mode. If the camera is disconnected from DC power, the battery will supply power for up to 2 minutes. If the camera is not reconnected to DC power within 2 minutes, the battery automatically deactivates and turns off the camera.

The battery operation mode does **not** automatically change to Capture when the camera is placed in Capture.

Capture Mode (Selectable): This mode allows the camera to operate in Capture mode for a short period without DC power. In the event of DC power loss, the battery will supply power to the camera either for 3, 5 or 10 minutes (Selectable). If the camera triggers within this time span, the battery operation mode automatically switches to Trigger Mode to protect the data. (Pertains to Stand- alone C321 only! The Miro JBox automatically triggers a C321J or C321 in

Capture mode in the event of power loss. This battery operation mode is disabled in the C321J).

Trigger mode (Default mode, always selected): This mode protects critical cines and ensures they are saved in the event of DC power loss. The battery activates when the camera is triggered. In the event of DC power loss, the battery will supply power to the camera until any of the following conditions are met:

- All auto save operations of Cines in RAM memory have been completed.
- All Cines in RAM memory have been deleted.
- The reset button on the back of the camera has been pressed.
- The battery voltage has reached the full discharge limit.

The camera then automatically turns off.

CHARGING

If the battery is not fully charged (at least 8.2 volts), it will automatically charge when the camera has DC power and is fully booted up. While charging, the camera's power draw is 22W, and the Battery LED on the back of the camera is Blue (if not armed) or Cyan (if armed).

The battery will not charge or stop charging under the following conditions:

- A battery fault is detected. In this case, the LED is Red.
- The battery will also stop charging automatically once it is fully charged. The LED light is off for a fully charged battery that is not armed.

A fully depleted battery requires 2.5 to 3 hours to fully charge.



Miro C321J Rear View



Miro C321 Rear View

The latest version of Phantom PCC software can be downloaded from the Vision Research website: www.phantomhighspeed.com/pcc

This manual covers the most commonly used functions. See the 'PCC Help' file for details of other settings.



PHANTOM CAMERA CONTROL (PCC) APPLICATION OVERVIEW

TOOLBAR

The '**Toolbar**' buttons provide quick access to the most frequently used functions. In the program, hover over a button to display a text box briefly describing its function.





The '**Help**' options provide valuable reference information, along with extensive documentation, relating to the software. Online tutorials can be found at *www.phantomhighspeed.com/tutorials.*

CONTROL TABS

PCC provides three control tabs: 'Live,' 'Play' and 'Manager.'

When first started, the '**Manager**' tab is selected. Connected cameras are selected for use and naming in this tab. It is also used to manage saved Cine files.

All camera control and capture parameters (sample rate, exposure time, etc.) are performed in the '**Live**' tab.

Reviewing, editing and saving of Cine files, either from the camera's internal RAM memory, installed Phantom CineMag V or external hard drive, are performed in the '**Play**' tab.

Live	Play	Manager	
0	◆s ×		
Car	eras 10277		
- Ö	12111		
B-C Flee	Contrast.		
P-	auto track		
	color matrix		
	New Upon		
10	neasurements		
	mage seach		
	tone curve		
- 22	chips		
- 2	justin blink		
- 22	Spay Bottle		
- 2	car2		
2	vapor trail ignit	on 29lps edited	
	vapor trail ignit	on 3200 ps edited	
	pvcpipe1_colo	r	

PHANTOM VIDEO PLAYER (PVP) APPLICATION OVERVIEW

PVP can be launched directly from the desktop shortcut or by the '**Video Out**' toolbar button in PCC.

PVP controls the video outputs connected to a video monitor or viewfinder only.

PVP provides the ability to view, capture, review, edit and/or save a Cine recorded into the camera's RAM, to a hard drive or installed Phantom CineMag. PVP is extremely effective when used with high-resolution cameras since most computers will not play captured raw files smoothly.

> PVP							-	0	×
Camera:	21111(211	11)		One	1	- 2	+5	Settings	
			Speed				***	Preference	
_	-	-	PlayBa	ck Speed		59	94	Help	
			Pa/Ne ⊠ Pi	c ly Each in	age 6.5	4(1/16		About.	
			Play M	ode				Spe	ed
		<u>"</u> U		nt To Ran	90 F	tay Al RA	м	60	T.
				ng Pong			1		
-	101	1997	Rash				-		
K	×	0.01	1 Free:	20000-06		Grase Al.		1 -	
-25477				8098					
-3321					1	1			
Card		Trigger			Save	to Resh		Save to file.	

The camera's video system and display settings are also set through PVP. The best video system for the project will vary based on the country you are in, what kind of video monitor is used and the required video resolution.

All available video systems for the connected cameras can be found in the '**Settings**' menu of PVP, along with production area and other video overlay controls.

CAMERA CONTROL VIA PCC

PCC provides the ability to select units for specific camera parameters by clicking the '**Preferences**' button at the bottom of the '**Manager**' tab.

Units can be set to commonly used values ('Presets') or can be customized using the pull-down selection lists. First time users should use one of the three '**Presets**.'

The '**Exp**' unit is probably the most important unit to be set. It specifies what unit to use when setting the exposure time. The other units to set are EDR (Extreme Dynamic Range, not compatible on all models) and PTF (Post Trigger Frames) covered later in this section.



SELECTING A CAMERA

Select the camera(s) to be controlled listed in the '**Manager**' tab, or select the camera(s) from the '**Camera**' pull-down list in the '**Live**' tab.

Once a camera is selected, a '**Preview**' panel will display to the left of the control tabs showing the current image being captured by the camera.

IMAGE PROCESSING

'Image Tools' provides extensive control over the look of the image, from color and contrast settings to image orientation and crop settings. The menu is accessed by clicking on the **'Image Tools**' toolbar button (the one that looks like an artist's palette).

The top of the 'Image Tools' window displays a 'Histogram.' This is a graphic representation of the pixel brightness levels of the displayed image. The left represents black, the right represents white and the height represents the proportionate number of pixels at that particular value. Unlike a waveform, the histogram's shape is not representative of the content—it is simply an averaging of the brightness values.

Below the histogram are controls which change image settings of the live images, recorded images and the video output.

Some of the variables include; brightness, gain, gamma, saturation, toe, white balance adjustments (Temp (K) and Tint), individual red, green and blue pedestal, gain and gamma values and tone curves.

At the bottom of the window is a '**Default**' button that restores all parameters except white balance, tone and color matrix to their default values.

The '**Default White Balance**' button restores white balance to the defaults on color cameras.



The Tone '**Reset**' button restores the image tone to the default values, and the Color Matrix '**Restore**' button returns the color matrix values to their default values.

Changes made here only affect the metadata of the Cine Raw file. They are applied in software but not "baked in."

Image Processing settings are no longer editable after saving to a format other than Cine Raw, so it is important to ensure they are set to values that produce the best looking image.

IMAGE DISPLAY

The 'Zoom Actual Size' toolbar button resizes the images being displayed in the 'Preview/ Playback' panel to their actual size.

The '**Zoom Fit**' toolbar button resizes the images to fit the panel. Images can also be zoomed to a specific magnification ratio by selecting a number from the pull-down list to the right of the '**Zoom Fit**' button.

WHITE BALANCE (COLOR CAMERAS)

Performing a 'White Balance' should be the first step in color adjustment. White Balance not applicable to monochrome cameras.

Right click on an area that resembles white or neutral area in the image, then click on the 'White Balance' pop-up window. It is not necessary to fill the frame with white—a small target can be used.



It is recommended to perform the White Balance after a CSR and on a white or gray object that is not fully saturated.

CAPTURE SETUP: CAMERA & CINE SETTINGS

Set Time: Synchronizes the timestamps embedded in the recorded image data to the computer's clock.

Bit Depth: Most Phantom cameras operate in 12-bit mode only.

Mode: Select the desired mode, binning or standard mode.

Partitions: Select the number of desired partitions (evenly divided memory segments) from the 'Partitions' pull-down menu. For basic camera setups, this should be set to one.

Lens Control: Available for Canon EF lenses only, for control of aperture and focus.

Backup & Restore: Allows for user and factory settings to be saved and recalled from the camera's memory.

Resolution: Sets the camera's acquisition resolution. There are several options in the pull-down menu. Alternatively, type in a value and the closest valid resolution will be set.

Sample Rate: Sets the acquisition frame rate in framesper-second (fps).

Exposure Time (shutter): Sets the exposure time in degrees, microseconds or percentage (this depends on how the PCC preferences are set).

Exposure Index: Sets the exposure index (Effective ISO) of the image by loading preset tone curves. Adjusting gamma, gain and other settings will contribute to the overall EI value, and this combined value is what should be used to determine lighting.

CSR (Current Session Reference): Closes the camera's internal shutter and resets the black point of every pixel for optimal image quality.

Image Range and Trigger Position: The slider represents the memory buffer, with the '**Duration**' indicated in seconds and the total number of frames available.

The trigger position is indicated in the **'Last**' pull-down menu and as the **'T**' slider along the timeline. The trigger position is exactly when the trigger signal will be detected in the Cine.

KEY ADVANCED SETTINGS

The "Start/End of Recording Actions" section provides options of actions that can automatically be performed at the start or end of each shot, including:

- Auto save to CineMag/Built-in Flash: This feature saves a user-specified portion of a clip to the internal Flash immediately after recording.
- Auto play Video Out: Begins playback after recording. The range marked under 'Auto play Video Out' affects both playback and saving to the internal Flash.

 Live
 Play
 Manager

 Camera:
 ■
 set to...

 12111
 ✓

 Image: Settings
 >

 Camera Settings

 Image: Setting

Continuous Recording Camera Info

Durrent Tile I	New 21 2017 17-54-02 Set
Time:	Time Time
6	Local O uto (GMT)
100 N/A	
ara. Inn	
Bit Depth 12	✓ Pattions 1 ✓
Sensor Acronit	tion Mode
HQ	✓ Set Mode
Lens Control	
Aperture	· Past+
~	Pocus 0
	Fast-
Use shin or Ch	for amaler focus steps.
Backup & resto	ire settings
Load	Save
V Cine Setting	
One 1	 ✓ m[™] set al.
-	
Name	
Name Description	2
Name Description	>
Name Description	2048 x 1952
Name Description	2048 x 1952 v
Name Description Resolution [Sample Rate [2048 x 1952 V 1000 V pps
Name Description Resolution Sample Rate [Exposure Time	2048 x 1952 V 1000 V pps 900 V µs
Name Description Resolution 2 Sample Rate 2 Exposure Time EDR	2048 x 1952 ~ 1000 ~ pps 1000 ~ µs 100 ~ µs
Name Description Resolution Sample Rate Exposure Time EDR	2048 x 1952 v 1000 v pps 2000 v µs 0 v µs
Name Description Resolution Sample Rate Exposure Time EDR Exposure Index	2048 x 1952 v 1000 v pps 2000 v µs 0 v µs Custom v
Name Description Resolution [Sample Rate [Exposure Time EDR Exposure Index EI + Image Proc	2048 x 1952 v 1000 v pos 500 v µa 0 v µa Custom v cessina 2500
Name Description Resolution 2 Sample Rate 2 Exposure Time EDR Exposure Index EI + Image Proc CSR	2048 x 1952 v 1000 v pps 1000 v µ8 0 v µ8 0 v µ8 Custom v tessina 2500 Low Light
Name Description Resolution Sample Rate Exposure Time EDR Exposure Index EI + Image Proc CSR Image Range a	2048 x 1952 v 1000 v pps 2000 v µs 0 v µs Custom v cessina 2500 Low Light and Trigger Position
Name Description	2048 x 1952 v 1000 v pps 500 v us 0 v us Custom v custom v Low Light and Trigger Position Lett: v
Name Description Resolution Exposure Time EDR Exposure Index EI + Image Proc CSR Image Range a -575 T	2048 x 1952 v 1000 v pps 900 v us 0 v us 0 v us Custom v costina 2500 Low Light Low Light Levt: v

• **Restart Recording:** When enabled, automatically restarts the recording process after the 'Auto' actions have been performed.

'External Sync' instructs the camera to use one of the following options:

- Internal: Camera uses its internal crystal oscillator to drive the frame rate.
- **External:** Camera uses externally supplied frame sync clock pulse to drive the frame rate. This can be used to synchronize two cameras together via F-Sync.
- IRIG: Camera uses an IRIG-B signal is supplied to drive the frame rate.
- Lock to Video: Frame rate is driven by the camera's current video rate. Fps will jump to the closest multiple of the current video rate (23.98, 24, 25, 29.97 or 30).
- **Sync to Trigger:** Instructs the camera to adjust its frame clock, upon detection of a trigger signal, to ensure all post trigger frames occur at the same moment in time from trigger when repetitive tests are required.

When using External Sync it is important to limit the top frame rate by ~1% to ensure the sync period is larger than 1/max_frame_rate + 0.1% or 200ns (whichever is larger), to account for frequency tolerance and jitter.

FLASH MEMORY

This displays the amount of free space and size (in Gigabytes) of the internal flash.

		Erace
1940.9	GB free of 2093.0 GB	
Dre	t recording to CineMag	

RECORDING A CINE

To begin recording to the camera's RAM click the red '**Capture**' button.

The red 'Capture' button changes to 'Abort Recording' and the green 'Trigger' button



is enabled when the camera is recording. The 'Abort Recording' button instructs the camera to stop recording, leaving the camera's RAM empty.

TRIGGERING THE CAMERA

Selecting the '**Trigger**' button instructs the camera to immediately stop recording when the '**Trigger Position**' is set to zero. If a value greater than zero is set, the camera will continue to record '**post-trigger**' frames until the user-specified value is met.

Using the camera's external trigger signal provides a more accurate trigger to the camera. This is also referred to as a 'hardware trigger.'

REVIEWING A CINE

Once the camera has completed recording a Cine in the camera's RAM or CineMag, it can be reviewed by selecting it from the '**Cine**' pull-down selection list in thePCC '**Play**' tab.

A previously saved Cine stored on the computer's hard drive can be opened using the '**Open File**' is toolbar button. (When used, it also places the file under the '**Cines**' group folder in the '**Manager'** tab.) Uve Play Manager
Crie:

I2111 > One 1

Files
viapor trail igntion 3200fps edited
procipie 1_color
Comerciae

I2111 > One F1
12111 > One F1
12111 > One F2
12111 > One F4

Use the '**Video Control**' buttons to review the Cine:





PERFORMING A QUICK SEARCH THROUGH A CINE

Quickly find the points of interest:

'Scroll' (scrub) through the clip using the slider or click anywhere on the timeline to jump to that point quickly.

'Jump' to the trigger frame by clicking on the **'T**' button, or jump to specific frames by entering the frame number into the jump **'#**' data entry field, then hit the enter key.

'Image Search.' The goal is to search or find an image change in the recording, based on the differences between image content. Right-click on the 'Play' button to begin the image search. Besides image content changes, 'Image Search' can also look for images that are tagged as 'Event' images.

-15725	-6803			15725
-15726	1 1 1	1.1	4	16726
Jump	H H	To #		
♥ Image Search Methods	1			
Difference >	Threshold:	2 %		
Changed areas:	10.00 ~ 3			
Search Step:	10			
Incremental s	earch			

EDITING AND SAVING A CINE

Using the following 'Video Control' buttons, locate the first image of the Cine to be saved and click the 'Mark-In' I button. Locate the last image of the Cine to be saved and click the 'Mark-Out' 1 button.

Click 'Play, Speed & Option' and enable (check) 'Limit to Range.'

Under the 'Video Control' buttons, click the play button and review the edited Cine.

Click the '**Save Cine...**' button to save the edited Cine to the computer's hard drive.

If you wish to save the clip to an internal Flash, click the down arrow to the right of the 'Save Cine...' button and select 'Save RAM Cine to Flash.'



WORKING WITH CINE FILES

The images recorded on the camera's RAM or Phantom CineMag are stored in a Vision Research proprietary RAW (uncompressed) file structure called a '**Cine**' file.

These Cine files can be converted to industry standard formats (ProRes, H264, DPX, DNG, TIFF, JPEG and more) with PCC software provided by Vision Research.

Phantom PCC and PVP software are compatible with Windows operating systems. There are third party solutions available for working with Phantom cameras in Mac OS X.

COMPATIBILITY WITH VIDEO EDITING PROGRAMS

Several popular video editing programs, such as DaVinci Resolve, have incorporated the Phantom Cine Raw file format into their software. This means Cine files do not have to be converted and no additional software is required.

Please test the footage with the program you choose before committing, as updates to the program or Cine file format can sometimes break compatibility. For this reason, it is important to know how to properly convert Cine Raw files using PCC.

CONVERTING CINE RAW FILES

Single Cine files can be converted by selecting the desired format from the '**Save as Type**' selection list in the '**Save Cine**' dialogue window.

The file formats above the separator line in the 'Save as Type' selection list are 'movie-

Re-saving a clip in the Cine Raw format can be useful for creating sub-clips with no loss in image quality or metadata.

like' formats (meaning the entire clip will be saved as a single file), while the formats below the line are image formats (meaning each frame of Cine will be saved as a sequence of images).

To convert a Cine to a '**movie-like**' format, select the desired format from the list, navigate to the destination folder, assign a file name to the clip and save.

Some valuable parameters can be found in the 'Advanced Settings' window, such as the particular codec.

Other formats, like .avi and .mp4, allow the compression ratio to be entered. The lowest compression is the default.

CONVERTING TO A SERIES OR STACK OF IMAGES

Beginning with PCC software version 3.5, the autoname functionality eliminates the need to add a special character to create a sequence of images. Use one of the presets in the save dialogue. Vision Research also recommends setting a project name, in the 'Auto-name settings' menu found in the top bar pull-down arrow next to the snapshot functions.

Cine Raw, *.cine	\sim
Cine, *.cine Cine JPEG, *.cci	^
Cine Raw * cine	
AVI. * avi	
Multipage TIFF, *tif	
H.264. *.mp4	
ProBes, * mov	
QuickTime Uncompressed, *.mov	
Windows BMP 24 images, *.bmp	
Windows BMP 8 images, *, bmp	
Windows BMP 4 images, *.bmp	
OS/2 BMP 4 images, * bmp	
OS/2 BMP 8 images, *.bmp	
OS/2 BMP 24 images, *.bmp	
PCX 1 images. *.pcx	
PCX 8 images. *.pcx	
PCX 24 images, *.pcx	
TGA 8 images, *tga	
TGA 16 images, *.tga	
TGA 24 images, *.tga	
TGA 32 images, *.tga	
TIFF 1 images, *.tif	
TIFF 8,24 images, *.tif	
TIFF 12,36 images, *.tif	
TIFF 16,48 images, *.tif	
LEAD images, *.cmp	
LEAD JFIF 4:1:1 images, *,iff	
LEAD JFIF 4:2:2 images, *.jff	
LEAD JTIF 4:1:1 images, *.jtf	
LEAD JTIF 4:2:2 images, *.jtf	
JPEG images, *.jpg	
JTIF images, *.jpg	
RAW images, *.raw	
DNG images, *.dng	
DPX images, *.dpx	
EXR images, *.exr	~

BATCH CONVERT

The 'Batch Convert Files' it toolbar button can be used to convert a single saved Cine file, or multiple saved Cine files, into any one of the supported file formats.

Use the shift and/or control keys to select the Cine files you wish to convert in the '**Open Cine**' dialogue window, then click the '**Open' button**.

Navigate to the destination folder in the '**Multifile Convert Destination**' dialogue window and select the file format.

PCC creates a separate folder for each Cine file, assigns the original file name and appends the appropriate image number and extension to each image.

Once the '**Convert**' button is clicked, a progress window appears for the duration of the conversion process. Each converted Cine will be placed in its own folder named after the original Cine file.

File: justin blink	20 %
Total 2/2	60%
	Close



PROGRAMMABLE I/O

Vision Research, Inc. | Miro C321/C321J/N5/JBox 2.0 Camera Manual



INTRODUCTION TO SIGNALS

Phantom cameras incorporate multiple hardware signals to interface to external devices.

Ports on Phantom Miro C321 cameras are programmable. The equipment to access these ports – the MiniBoB, Control BoB, and Junction Box – have ports that are either fixed or programmable. The Programmable I/O ports have a default setting which will be set after a factory reset or by using the 'signal default' function in PCC.

Port Map for Programmable I/O Access

MINIBoB								
Port Label	Default setting	Programmable I/O Capability	Programmable IO Port in PCC					
Trigger	Trigger	Fixed	Port 1					
TC In	TC In	Fixed	Port 2					
TC Out	TC Out	Prog. I/O	Port 3					
Aux 1	Strobe	Prog. I/O	Port 4					
Aux 2	Ready	Prog. I/O	Port 5					
	CONTR	OL BoB						
Port Label	Default setting	Programmable I/O Capability	Programmable IO Port in PCC					
Trigger	Fixed	Fixed	Port 1					
TC In	Fixed	Fixed	Port 2					
Ready	Ready	Prog. I/O	Port 5					
TC Out	TC Out	Prog. I/O	Port 3					
Aux Out	Strobe	Prog. I/O (out)	Port 4					
Aux In	N/A	Prog. I/O (in)	Port 4					

Programmable I/O incorporates a pulse processor. In addition to assigning different signals, the signal characteristics can be modified to better interface with external devices. In most cases, the signal polarity, filter time, delay and pulse width and edge (rising versus falling) can be set. Setting these characteristics is referred to as 'Pulse Processor Control.'

PCC INTERFACE

The 'Camera Signals' menu, located in the 'Live' tab, provides access to and control over these Programmable I/O signals.

V Camera Signals 1 Trigger ~ 0 Faling Edge Filter: 32 µa Delay: 0 µa 2 Time Code In Fixed Port Programmable I/O: Default All P3 Frame Sync I/O ~ 0 Filter: 0 µ8 P5 Ready Width: 0 µs Delay: 0 µs Omba P4 Strobe Frame Sync I/O Time Code Fiber: 0 µs Multi-Strobe Width: 0 µs Delay: 0 µs Autotrigge Software Trigger Recording P5 Ready ~ 0 Filter: 0 up A.e. Event In Memory Gate In Width: 0 µs Delay: 0 µs Frat Pretrigger In Ima Runstop In Cor Austrager In Can Core Event Core Memory Gate P6 Time Code ~ Q Fiber: 0 µ8 Width: 0 µs Delay: 0 µs Core Frame Sync Core Pretrigger Core Runstop Core Austrigger te Trioce

A summary of the current signal settings can be found under the pulldown menu associated with each port. By clicking '**Default All**,' PCC will reassign the factory-assigned defaults for the Programmable I/O signals.

PULSE PROCESSOR CONTROL

After selecting the signal assignment, further configuration is possible by clicking the gear symbol next to each port. The '**Pulse Control**' menu is opened, as shown on the next page.

A graphic representation of the signal behavior is also displayed. However, this is **not to scale** and should be used as reference only. Use of this feature requires an oscilloscope to truly visualize the signals and the subsequent changes with each adjustment.

Programmable I/O	Port 4 Pulse Con	trol		×	Programmable I/O P	Port 4 Pulse	Control		×
Signal Name: Fra Invest Wodth Delay Filter	e Sync I/O Faling Edge 2.000 1.000	2 2	Unte Orme ⊛µa Orme O Occka		Signal Name: Fran	e Sync I/O Falling E 0.000 0.000	dor M M M	Unite ○ mi ● µi ○ ni ○ Ciocka	
Frame Sync I/O (Processed)	-> Deay	4- With A-	Shiph Not To Scale	-	Frame Sync I/O (Processed) Default	-		Orașh Nor To Soare	ncel

Input: When a signal is generated by the camera, it serves as an input to the 'pulse processor.' In this context, the term 'input' does not represent an externally generated signal.

SUMMARY OF PULSE PROCESSOR SETTINGS

Invert: Inverts the signal at the output of the pulse processor.

Falling: Selects 'Falling Edge' mode for the pulse processor. This mode is only relevant if the 'Width' is also specified. When the 'Falling' token is present together with 'Width,' the pulse processor will generate a negative pulse, triggered from the negative edge of the input signal.

Width (Pulse Width): When a 'Width' token is present, a defined-length pulse is generated, which starts after the specified 'Delay' and after the active edge of the 'input' signal. The length of the pulse is specified in microseconds (as a floating point number) and internally converted to pixel clock units. The maximum pulse width is at least 10 seconds. However, if the period of the 'input' signal is lower than the selected width, the latter is dynamically clamped to the signal period. The minimum pulse width is one pixel clock.

Delay: Delays the output pulse by the specified time in ms, µs or camera clock multiples. If the 'Width' token is not present, both edges of the signal are delayed by the same amount. If present, the delay is measured from the rising edge of the input* signal unless the 'Falling' token is present, in which case the delay is measured from the falling edge of the input. The delay time is specified in microseconds and is internally converted to pixel clock units.

Filter (Filter Time): When a 'Filter' token is present, the 'input' of the pulse processor is filtered through an edge filter of the specified time. The time of the filter can be between 0 and 1 second. In order for the output of the filter to be asserted, the 'input' signal must be continuously deasserted for the same duration. The edges of the 'input' are thus delayed by the specified filter time (for a 'clean' input pulse). Filtering is applied before and independently of the delay and duration. The filter time is specified in microseconds and is internally routed to pixel clock units.

Pixel Clock: The period of the pixel clock is the basic time interval for all camera timing.

Please note that pulse processors can sometimes generate pulses too short for the output drivers to switch and the processed signal should be verified with an oscilloscope before use.

PROGRAMMABLE SIGNALS & DESCRIPTIONS

The settings as described are the signal's default state prior to processing.

Ready: An isolated open collector output with 1k pull-up signal (active high). 'Ready' is asserted when the camera goes into 'Capture' mode and is de-asserted either when the Cine is triggered or when recording is completed. 'Ready' changes synchronously with frame capture (at the end of each exposure), so in 'External Sync' mode it will not change until F-Sync pulses are received.

Strobe: An isolated open collector output signal with 1k pull-up. When asserted (low) 'Strobe' indicates that the electronic shutter is open. Strobe is low for the duration of the exposure.

F-Sync: The only signal can be set as an output or input. By default, it is output (sync-internal). Output signal is a frame sync pulse from the camera's frame rate generator. A short (few

hundred ns depending on camera model) negative pulse with the falling edge is used as timing reference. Input signal is active on falling edge. (Default state is high.)

TC-Out: A positive polarity time code signal. Normally an unmodulated (dc-shifted) IRIG-B (at RS-232 levels) which follows the internal time base of the camera. It is recommended not to process the 'TC-Out,' since a processed signal will not represent an accurate time code.

Auto-Trigger: Used to output a hardware trigger signal or pulse with the duration.

Software Trigger: An active high-output signal (pulse) generated as a result of the trigger protocol command.

Recording: An active high-output signal indicating active recording into a RAM partition.

Event In: If the input is sampled low at the end of an exposure, an 'E' (Event) bit in the frame's timestamp is set.

Memory Gate In: If the input is sampled low at the end of an exposure, the corresponding frame is skipped from storage to RAM.

Pre-trigger: An active low-input (default high) signal. Keeping this signal low for enough time (10–500 ms or until 'Ready' signal goes high) will make the camera start recording, if it has an available RAM partition.

Aux Trigger: An input signal active on the rising edge (default high). An alternative trigger input that can be processed through the pulse processor and assigned to different ports.

THE CORE SIGNALS

Core signals are copies of externally generated signals, routed through the camera and output to assigned ports. Unless indicated otherwise, feedback is taken after any pulse processor for the input. Core signals can be pulse-processed before being output.

Core Event: Feedback output from the 'Event In' signal.

Core Memory Gate: Feedback output from the 'Memory Gate In' signal.

Core Frame Sync: Feedback output from the 'F-Sync In' signal. The feedback is taken after any pulse processor for the output, but before the delay element.

Core Pre-trigger: Feedback output from the 'Pre-trigger' signal.

Core Aux Trigger: Feedback output from the 'Aux Trigger In' signal.

Core Trigger: Feedback output from the main 'Trigger' input. The feedback is taken before the trigger signal is affected by the trigger polarity, filter or delay settings. 'Core Trigger' can be used like the 'Trigger Out' signal.





Phantom Miro Junction (JBox) Box

MIRO JUNCTION BOX 2.0 (JBOX 2.0)

The Miro Junction Box 2.0 (JBox 2.0) is modular and flexible: with six camera ports, it can host up six cameras or dedicate any port to uplink to another JBox in a daisy chain or branch to expand the number of cameras in a configuration. It is compatible with all Miro C Series, N5, and VEOs, as well as Miro M, R, and LC Series.



Every JBox requires 20 - 36 VDC of power to power the cameras connected to it. Cameras connected to a JBox receive power through the JBox and should not be connected to an additional power source. The number of cameras and type of cameras connected to a single JBox determines how much wattage and voltage is required.

JBOX 2.0 POWER Max cable length of each camera = 30M

Option	Port 1	Port 2	Port 3	Port 4	Port 5	Port 6	Power Req.
1	Miro C	280W					
2	Miro R	Miro C	280W				
3	Miro R	Miro R	Miro C	Miro C	Miro C	Miro C	280W
4	Miro R	Miro R	Miro R	Miro C	Miro C	Miro C	280W
5	Miro R	Miro R	Miro R	Miro R	Miro C	Miro C	280W
6	Miro R		280W				
7	Miro R	320W, 32V					
8	VEO S	Miro C	320W, 32V				
9	VEO S	VEO S	Miro C	Miro C	Miro C	Miro C	320W, 32V
10	VEO S	VEO S	VEO 4K	Miro C	Miro C	Miro C	320W, 32V
11	VEO S	VEO S	VEO S	VEO S			320W, 32V
12	VEO S	VEO S	VEO 4K	VEO 4K			320W, 32V
13	VEO S	Miro R	480W				
14	VEO S	VEO S	Miro R	Miro R	Miro R	Miro R	480W
15				Miro R	Miro R	Miro R	480W
16	VEO S	VEO S	VEO 4K	VEO 4K	Miro R	Miro R	480W
17	VEO S	VEO S	VEO 4K	VEO 4K	Miro C	Miro C	480W



CAMERA CONTROL

The JBox offers 2 methods to control the cameras:

- 1. Via Gb Ethernet, controlled directly by a computer with PCC or third-party software.
- 2. Via the System Port and a system cable to access Trigger, IRIG, Strobe, FSYNC, Ready control, and Ethernet. The cable may be spliced or connected to a Control Break-out Box (Control BoB). More detail about the Control BoB is in the Accessories Section.

SIGNAL MANAGEMENT

PCC: Aux 2 should be set to "Ready" for all cameras; Aux 3 should be set to "IRIG" for all cameras.

Ethernet: The JBox has an RJ45 Ethernet connection

Trigger: Does not require a specific set-up in PCC. The JBox accepts a Trigger signal either through the System cable, or through the Trigger BNC.

Synchronization by IRIG: Set all cameras "Sync to IRIG" in PCC.

The JBox accepts an IRIG signal from an external source through the System cable. If no external source is present, the JBox will select one of the attached cameras as the IRIG source.

Synchronization by FSYNC: In PCC, Aux 1 should be set to "FSYNC" for all cameras.

Set all cameras except one to "Sync to External" in PCC. Set the one remaining camera to "Sync to Internal". This one camera may be in any port.

If using a camera FSYNC do not use any external FSYNC signals.

The JBox accepts an FSYNC signal either via the system cable, or via the Aux 1 BNC, or via any camera connected to the JBox.

AUXILIARY FUNCTIONS

Strobe: In PCC, Aux 1 should be set to "Strobe" for all cameras.

The JBox automatically selects one of the cameras to emit a strobe signal. The signal is output to both the Aux BNC and the System cable. The cameras connected to the JBox capture images synchronously based on the IRIG setting. Set the same frame rate and exposure for all cameras, to ensure they all capture images simultaneously.

It is not possible to sync via FSYNC and use Strobe at the same time. **Event:** In PCC, Aux 1 should be set to "Event/Memgate" for all cameras.

The Event signal may be input into either the Aux BNC or the System cable.

Ready: Aux 2 in PCC should still be set to "Ready". All cameras connected to the JBox

automatically emit a Ready signal when in capture mode.

All cameras emit a Ready signal. If all cameras are Ready, a Ready signal is output to the Ready BNC and the System cable. In a Daisy Chain configuration of connected JBoxes, the Ready signal flows from the slave JBoxes to the master JBox (closest to the control room). This is the opposite direction from the other signals. Only the Ready signal from the master JBox represents the entire system.

ADVANCED JUNCTION BOX 2.0 CONFIGURATIONS

Multiple Junction Boxes may be connected via system cables to create "daisy chain" or "tree branch" configurations. A camera port of the first JBox is connected via system cable to the System port of the next JBox, and so on to build either a chain, branch or combination configuration.

Power: In a multiple JBox configuration,⁴ each JBox requires a power source to power the camera connected to it.



Signal Flow: Signals begin in the Master JBox (Closest to the control room), then flow to the other JBoxes in the configuration.

- Ethernet: Ethernet should be connected to just the master JBox. Ethernet will flow to the other JBoxes.
- Trigger: Trigger should be connected to the master JBox. It will flow to the other JBoxes.
- IRIG: The IRIG signal will be generated from the first camera in the master JBox, and flow to the other JBoxes.
- FSYNC, Event, MemGate Inputs: These inputs should be connected to the master JBox, either through the AUX BNC or the System cable. They will flow to the other JBoxes.
- FSYNC, Strobe Outputs: These outputs will be generated by the first camera in the master JBox. They will flow to the other JBoxes.
- Ready: The ready signal is the only signal that flows from the farthest JBox back to the master JBox. A JBox that is not the master may indicate that its cameras are ready; however only the Ready signal from the master JBox represents the entire system.



CONTROL BREAKOUT BOX

The Control Breakout Box (Control BoB) offers a sleek and simple way to connect the required signaling to a Junction Box configuration without splicing cables. With ports for IRIG, Ethernet, Ready, Strobe, and FSYNC, the Control BoB is a clean and convenient solution to manage signaling. It measures 7.4 x 2.4 inches (187mm x 60 mm).

It functions using 7-32VDC, and has two power ports. Choose either the rugged Amphenol port with a customer supplied power supply for use in rugged environments, or the barrel port with a supplied power supply for non-rugged use. Each Control BoB must have a power supply and draws 0.15Amps. The Control BoB cannot power the Junction Box or a camera.



MIRO CONTROL BoB

A TRIGGER

Trigger (input): When a TTL pulse (rising/falling edge) is detected, the camera triggers.



Ready: Outputs a TTL signal (high) to indicate that all cameras connected to the Miro JBox 2.0 are in Capture Mode. Signal can be set to go low at trigger or at the end of recording.



TC In (input): Connects to an IRIG receiver to accept an unmodulated signal to drive the Miro Junction Box IRIG signal generation.



Aux IN: The following inputs can be used in this port:

- EVENT (input): When the Event signal is active, frames are tagged with an Event marker (as metadata). These events can be searched or referenced during playback.
- MEMGATE (input): When Memgate signal is active the camera stops recording into its internal memory (frames are discarded).
- F-SYNC (input): Connect an external source, including the F-Sync from a second Phantom camera, to drive the camera's frame rate. Use in combination with Sync: External in the External Sync menu.



TC Out (output): Connect to a second Phantom camera, to drive the camera's frame rate. Use in combination with Sync: Lock to IRIG in the External Sync menu.

AUX OUT

Aux OUT: The following outputs can be used in this port:

- STROBE (output): Signal goes low for the duration of each frame's exposure.
- F-SYNC (output): FSYNC output from active camera connected to the Miro Junction Box.

G POWER 7-32VDC

Power Connector: Amphenol connector for rugged applications for customer supplied power supply 7-32VDC, 500mA.



Ethernet: Ruggedized RJ45 connector connects GigE to a control unit computer / laptop for camera control communication.



Power Connector: Barrel connector for power supply 7-32VCD, 500mA.



System Connector: Same Fischer connector as System port on JBox 2.0. The Control BoB must connect to the System Port (not a Camera Port) of the JBox 2.0 to function as designed.

CABLES FOR JUNCTION BOX 2.0



System cable, Adapter cable, Short cable to link JBoxes

System Cables: Phantom system cables have differential signals for signal integrity over longer distances. They are capable of transporting: Power, Ethernet, and IRIG, FSYNC, Strobe, Ready, Memgate and Event signals. They have a rugged casing and the same Fischer connect at each end, one in a 90° angle and the other in a straight format. Either end connects directly to any port on the JBox 2.0. The other end connects directly to either:

- Miro C321J, N5 N-JB Base, C320J, or C210J
- Adapter cable

Adapter Cables: Phantom adapter cables connect VEOs and non-J Miro C cameras to a System cable for connection to a JBox 2.0. The adapter cable is approximately 0.6M (2Ft) and connects to the power and capture ports of the camera and mates with a system cable. The cameras it supports are:

- All VEO S-Models
- Miro C321, N5 N-B Base, C320, C210

JBox Connector Cable: This small cable is perfect for connecting 2 JBoxes that are immediately next to each other. It is a System cable that is only 0.6M (2Ft) long with both connectors at 90° angles. One connector attaches to a camera port of the master JBox, and the other to the System port of the next JBox.

MIRO MINI BREAK OUT BOX



The Miro C321 Capture port is compatible with a Phantom Miro Mini Break Out Box (Miro MiniBoB).

This connector is not used for the Miro C-Series.



TC In (input): Connects to an IRIG receiver to accept a modulated or unmodulated signal to drive the camera's frame rate.

C TC OUT

TC Out (output): Connect an a second Phantom camera, to drive the camera's frame rate. Use in combination with Sync: Lock to IRIG in the External Sync menu.

D AUX 1

Aux 1 (input / output switchable):

- STROBE (output): Signal goes low for the duration of each frame's exposure.
- EVENT (input): When the Event signal is active, frames are tagged with an Event marker (as metadata). These events can be searched or referenced during playback.
- MEMGATE (input): When Memgate signal is active the camera stops recording into it's internal memory (frames are discarded).
- F-SYNC (input / output): Connect an external source, including the F-Sync from a second Phantom camera, to drive the camera's frame rate. Use in combination with Sync: External in the External Sync menu.

Aux: 2: (output switchable):

- READY (output) When signal is high it indicates that the camera is in capture mode. Using PCC, signal can be set to go low at trigger or at the end of recording.
- STROBE (see Aux 1 description above).
- **TRIGGERTrigger (input):** When a TTL pulse (rising/falling edge) is detected,
the camera trigger isolated, \leftarrow 1V deassert, 3 to 24V assert.

E AUX 2 Au

MOUNTING ACCESSORIES



Camera / lens mount: A camera/lens mount makes mounting the camera easier and also provides stability and protection for the lens.









= 16.0 🗕

46.3 -

--- 7.5



MIRO C321 BOTTOM VIEW



MIRO C321 RIGHT / LEFT





MIRO C321 FRONT / BACK





N5 & BASES TOP / BOTTOM / SIDE























Use these schematics to build custom cables at your own risk. Miswired cables can cause serious damage to the camera, which is not covered under warranty. Vision Research recommends only using cables supplied by Vision Research.

These pin-out diagrams refer to the connector on the camera or Miro Junction Box body. Part numbers indicated are for the cable's connector.

MIRO C321 CAPTURE CONNECTOR

Capture port (for capture cable and MiniBoB) 12-pin Fischer Part # S-1031-Z012-130



PIN	NOMENCLATURE/FUNCTION
1	IOGND / Signal Ground
2	IOGND / Signal Ground
3	IOGND / Signal Ground
4	TRIGGER / Isolated Input. Active low. Can be activated by a switch to ground. The trigger pulse needs to be at least 3 microseconds long. Capture cable = Red
5	AUX2 (READY) / Isolated collector output with 1k pull-up. Capture cable = Green
6	AUX1 / Isolated collector output. Switchable between STROBE, EVENT, MEMGATE and F-SYNC. Capture cable = Black
7	TCIN / Timecode Input can accept IRIG-B and SMPTE standards. Capture cable = White
8	GND (reference for Pin 9)
9	Not used
10	POWER / Nominal power supply voltage is +24VDC (acceptable range is 16-36VDC)
11	POWER / (see Pin 10)
12	TCOUT / Timecode Output for IRIG-B or SMPTE timecode. IRIG Swings to RS-232 levels of +/- 9V. Capture cable = Blue

Vision Research, Inc. | Miro C321/C321J/N5/JBox 2.0 Camera Manual

CONNECTOR PIN-OU

MIRO C321 GIGE ETHERNET CONNECTOR

Ethernet Port

8-pin Fischer Part # SS 103 A058-130



PIN	NOMENCLATURE/FUNCTION
1	ETHRXP / 10/100/1000BASE-T Ethernet Receive (positive)
2	ETHRXN / 10/100/1000BASE-T Ethernet Receive (negative)
3	ETHTXP / 10/100/1000BASE-T Ethernet Transmit (positive)
4	ETHTXN / 10/100/1000BASE-T Ethernet Transmit (negative)
5	MDI2P /10/100/1000BASE-T Media Dependent Interface 2 (positive)
6	MDI2N / 10/100/1000BASE-T Media Dependent Interface 2 (negative)
7	MDI2P / 10/100/1000BASE-T Media Dependent Interface 3 (positive)
8	MDI3N / 100/1000BASE-T Media Dependent Interface 3 (negative)

MIRO C321 +16-32VDC CONNECTOR

Power Port

6-pin Fischer Part # SS 103 Z 056 130



PIN	NOMENCLATURE/FUNCTION
1	PGND / Power Ground
2	+VINBF / The +VDC In (Battery Freed) connector provides +24VDC (Direct Current) positive power to the Phantom camera.
3	RxD1 / RS-232 Receive Data 1
4	TxD1 / RS-232 Transmit Data 1
5	UNUSED
6	GND / Ground

MIRO C321J SYSTEM CONNECTOR

System Port (for Miro Junction Box)

27-pin Fischer Part # SS 105A 102 130



PIN	NOMENCLATURE/FUNCTION
1	GND / Power Ground
2	GND / Power Ground
3	GND / Power Ground
4	TRIGIN- / Trigger input (neg) from J-Box
5	READYOUT- / Ready / Strobe output (neg) to J-Box
6	IRIGIN- / IRIG input (negative) from J-Box
7	AUXIN+ / Strobe / Event / Memgate / Fsync input (positive) from J-Box
8	CBLDET / Cable detect. Connect to GND in connector
9	PWRIN_A / +20-32 VDC input from J-Box
10	MDI2- / Gigabit Ethernet
11	MDI1- / Gigabit Ethernet
12	MDI0- / Gigabit Ethernet
13	AUXOUT+
14	TRIGIN+ / Trigger input (pos) from J-Box
15	READYOUT+ / Ready output (pos) to J-Box
16	IRIGOUT+ / IRIG output (pos) to J-Box
17	IRIGOUT- / IRIG output (neg) to J-Bax
18	IRIGIN+ / IRIG input (pos) from J-Box
19	AUXIN- / Strobe / Event / Memgate / Fsync input (neg) from J-Box
20	AUXOUT-
21	PWRIN_B / +20-32 VDC input from J-Box
22	PWRIN_C / +20-32 VDC input from J-Box
23	MDI3+ / Gigabit Ethernet
24	MDI3- / Gigabit Ethernet
25	MDI2+ / Gigabit Ethernet
26	MDI1+ / Gigabit Ethernet
27	MDI0+ / Gigabit Ethernet

C321J REMOTE CONNECTOR

Remote Port (for Phantom RCU) 6-pin Fischer Part # SS 103 Z 056 130



PIN	NOMENCLATURE/FUNCTION
1	PGND / Power Ground
2	+VINBF / The +VDC In (Battery Freed) connector provides +24VDC (Direct Current) positive power to the Phantom camera.
3	RxD1 / RS-232 Receive Data 1
4	TxD1 / RS-232 Transmit Data 1
5	UNUSED
6	GND / Ground

MIRO JBOX 2.0 POWER CONNECTOR

Power Port 4-pin Amphenol Part # PT06A-12-4S



PIN	NOMENCLATURE/FUNCTION
А	+24VB / connect to a +24VDC battery running parallel to power source
В	CHGND / Chassis Ground
С	GND / Power Ground
D	+24VDC/ +24VDC (direct current) to J-Box runs parallel with an optional +24VDC battery

MIRO JBOX 2.0 ETHERNET CONNECTOR

Ethernet Port RJ-45



PIN	NOMENCLATURE/FUNCTION
1	MDI2+ / Gigabit Ethernet
2	MDI2- / Gigabit Ethernet
3	MDI3+ / Gigabit Ethernet
4	MDI3- / Gigabit Ethernet
5	MDI1+ / Gigabit Ethernet
6	MDI1- / Gigabit Ethernet
7	MDI0+ / Gigabit Ethernet
8	MDI0- / Gigabit Ethernet

MIRO JBOX 2.0 SYSTEM CONNECTOR

System Port (for Miro C321J)

27-pin Fischer Part # SS 105A 102 130



PIN	NOMENCLATURE/FUNCTION
1	GND_A / Signal ground for AUXIN and READYOUT
2	UNUSED
3	GND_B / Signal ground for IRIGIN and IRIGOUT
4	TRIGIN- / Trigger input (neg) to J-Box
5	READYOUT- / Ready / Strobe output (neg) from J-Box
6	IRIGIN- / IRIG input (negative) to J-Box
7	AUXIN+ / Strobe / Event / Memgate / Fsync input (positive) to J-Box
8	UNUSED
9	UNUSED
10	MDI2- / Gigabit Ethernet
11	MDI1- / Gigabit Ethernet
12	MDI0- / Gigabit Ethernet
13	AUXOUT+
14	TRIGIN+ / Trigger input (pos) to J-Box
15	READYOUT+ / Ready output (pos) from J-Box
16	IRIGOUT+ / IRIG output (pos) from J-Box
17	IRIGOUT- / IRIG output (neg) from J-Box
18	IRIGIN+ / IRIG input (pos) to J-Box
19	AUXIN- / Strobe / Event / Memgate / Fsync input (neg) to J-Box
20	AUXOUT-
21	UNUSED
22	UNUSED
23	MDI3+ / Gigabit Ethernet
24	MDI3- / Gigabit Ethernet
25	MDI2+ / Gigabit Ethernet
26	MDI1+ / Gigabit Ethernet
27	MDI0+ / Gigabit Ethernet

MIRO JBOX 2.0 CAMERA CONNECTOR

Camera Port (for Miro C321J)

27-pin Fischer Part # SS 105A 102 130



PIN	NOMENCLATURE/FUNCTION
1	GND / Power Ground
2	GND / Power Ground
3	GND / Power Ground
4	TRIGIN- / Trigger input (neg) to camera
5	READYOUT- / Ready / Strobe output (neg) from camera
6	IRIGIN- / IRIG input (negative) to camera
7	AUXIN+ / Strobe / Event / Memgate / Fsync input (positive) to camera
8	CBLDET / Cable detect. Connect to GND in connector.
9	PWRIN_A / +20-32 VDC input to camera
10	MDI2- / Gigabit Ethernet
11	MDI1- / Gigabit Ethernet
12	MDI0- / Gigabit Ethernet
13	AUXOUT+
14	TRIGIN+ / Trigger input (pos) to camera
15	READYOUT+ / Ready output (pos) from camera
16	IRIGOUT+ / IRIG output (pos) from camera
17	IRIGOUT- / IRIG output (neg) from camera
18	IRIGIN+ / IRIG input (pos) to camera
19	AUXIN- / Strobe / Event / Memgate / Fsync input (neg) to camera
20	AUXOUT-
21	PWRIN_B / +20-32 VDC input to camera
22	PWRIN_C / +20-32 VDC input to camera
23	MDI3+ / Gigabit Ethernet
24	MDI3- / Gigabit Ethernet
25	MDI2+ / Gigabit Ethernet
26	MDI1+ / Gigabit Ethernet
27	MDI0+ / Gigabit Ethernet

CONTROL BREAK OUT BOX (BoB) SYSTEM CONNECTOR

System port (for J-Box) 27-pin Fischer Part # SS-105A-102-130



PIN	NOMENCLATURE/FUNCTION
1	UNUSED
2	GND/ Ground
3	UNUSED
4	TRIGIN- / Trigger input (neg) to J-Box
5	READYOUT- / Ready / Strobe output (neg) from J-Box
6	IRIGIN- / IRIG input (negative) to J-Box
7	AUXIN+ / Strobe / Event / Memgate / Fsync input (positive) to J-Box
8	UNUSED
9	UNUSED
10	MDI2- / Gigabit Ethernet
11	MDI1- / Gigabit Ethernet
12	MDI0- / Gigabit Ethernet
13	AUXOUT+
14	TRIGIN+ / Trigger input (pos) to J-Box
15	READYOUT+ / Ready output (pos) from J-Box
16	IRIGOUT+ / IRIG output (pos) from J-Box
17	IRIGOUT- / IRIG output (neg) from J-Box
18	IRIGIN+ / IRIG input (pos) to J-Box
19	AUXIN- / Strobe / Event / Memgate / Fsync input (neg) to J-Box
20	AUXOUT-
21	UNUSED
22	UNUSED
23	MDI3+ / Gigabit Ethernet
24	MDI3- / Gigabit Ethernet
25	MDI2+ / Gigabit Ethernet
26	MDI1+ / Gigabit Ethernet
27	MDI0+ / Gigabit Ethernet

CONTROL BoB POWER CONNECTOR

4-pin Amphenol Part # 06A84S0



PIN	NOMENCLATURE/FUNCTION
А	UNUSED
D	+24VDC – Provide +24VDC positive power to Control BoB
С	PGND – Power Ground
В	UNUSED



Miro C321J Rear View



Miro C321 Rear View

LED INDICATORS

MIRO C321/C321J ETHERNET LED

A ETHERNET LED

COLOR	ETHERNET STATE
Green	Ethernet Link
Amber	Ethernet Activity

MIRO C321/C321J SYSTEM/STATUS LED

B SYSTEM / STATUS LED

COLOR	CAMERA STATE
White	Camera Booting
Green	Preview Mode
Red	Capture Mode

MIRO C321/C321J BATTERY LED

C BATTERY LED

COLOR	BATTERY STATE
Off	Battery fully charged, but not armed
Green	Battery armed
Blue	Battery charging or DC power removed
Red	Battery charging fault
Cyan	Battery armed and charging or DC power removed
Purple	Battery charging or DC power removed and charging Fault
Amber	Battery armed and charging fault
White	Battery armed, charging or DC power removed, and charging fault

MIRO JBOX 2.0 STATUS INDICATORS



Miro JBox 2.0

A POWER 20-36VDC INDICATOR

COLOR	POWER STATUS
Green	Power Good
Blue	Under voltage
Red	Overvoltage or Overcurrent

B READY INDICATOR

COLOR	READY STATUS
Green	Ready
Red	Not Ready

C TRIGGER INDICATOR

COLOR	TRIGGER STATUS
Green	BNC Trigger Asserted
Blue	Uplink / Expand Trigger Asserted

D AUX INDICATOR

COLOR	AUX STATUS
Green	Output
Blue	Input

E ETHERNET INDICATOR

COLOR	CONTROLLER ETHERNET STATUS
Green	EtherLink Active
Yellow	Ethernet Activity

F SYSTEM INDICATOR

COLOR	SYSTEM STATUS
Green	EtherLink Active
Yellow	Ethernet Activity

G CAMERA STATUS INDICATOR

COLOR	CAMERA STATUS
Green	Ready
Red	Not Ready

H CAMERA ETHERNET INDICATOR

COLOR	CAMERA ETHERNET STATUS
Green	EtherLink Active
Red	Ethernet Activity

MIRO C FAQS

What is the difference between the Miro C321Jand Miro C321?

The difference is in the connections. The Miro C321J must be operated through a system cable to the Miro Junction Box and does not have a separate power connector. It is the perfect, lowest cost solution for a system of multiple connected cameras. The Miro C321 has the same 3 connectors as other Miros. It can be powered and operated independently, or, with a Miro C Capture cable, it can also connect to the Miro Junction Box.

Otherwise, the Miro C321J and Miro C321 are the same.

Set Default CSR?

The Miro C's do not have a mechanical shutter. This allows them to be as rugged as they are but can make a CSR for good image quality inconvenient because you need to cover the lens. When you turn the camera off, the CSR returns to the factory default.

For customers that rarely change their camera settings, the "Set new CSR default" feature provides a convenient alternative to this process. You must operate the camera at full resolution, for example 1920 x 1080 for the C321. You may then set the CSR for the frame rate and exposure time you are using. That CSR will remain through turning the camera off and on, for as long as you don't change that setting. Always run at 1000 fps and a certain exposure, for example? Set CSR once, get great image quality, and don't worry about it again, no matter how many times you turn the cameras on and off.

Can I use my F-mount lens with the Miro C-Series cameras?

Yes, with a converter. The Miro C's come with a C lens mount that can be reversed to become a CS lens mount. To use an F-mount lens, you can purchase a converter (part number VRI-FMNT-CMNT).

Does the MiniBoB work with these cameras? What signals are available?

The Miro C321J does not use a MiniBoB, because it receives signaling from the JBox 2.0 through the System cable. The Miro C321 can use the MiniBoB.

The signals available to both cameras are:

Trigger F-Syn Memgate Ready IRIG In Strobe IRIG Out Event

What is the worst case power draw of the camera?

The cameras draw 22W while the batteries are charging.Otherwise, they draw only 12W during operation.

MIRO JBOX 2.0 FAQS

Which camera can I use with the Miro Junction Box?

The Miro Junction Box (JBox 2.0) accepts all Miro and VEO S Model cameras – The Miro C321J uses a System Cable, all other Miro and VEO S cameras require an Adapter Cable. Please note that the Miro JBox cannot accept other Phantom cameras.

If boxes are connected together, how many power sources are needed?

Each JBox 2.0 needs its own power source, such as a power supply, even if they are connected together.

BASIC TROUBLESHOOTING

There are fixed black spots on the image, particularly at small apertures.

This is most likely caused by dust particles stuck on the sensor or OLPF (optical low pass filter) surface. The best way to check is to remove the lens and look at the glass surface with a bright light source. Vision Research recommends not to use canned air or cotton swabs to clean the sensor surface. It is safer to use a manual bulb-style air blower and/or sensor cleaning brush for removing loose dust particles.

Additional tips for cleaning sensors can be found in the maintenance section of this manual.

TROUBLE SHOOTING

The live images look scrambled and the frame rate can't be set.

If the image is non-responsive and scrambled, the camera may be set to an external sync without a sync source connected. In the PCC \rightarrow Live \rightarrow Advanced Settings menu, check the External Sync setting to ensure that 'Internal' is selected.

Restoring the factory defaults.

If the camera is stuck in an unusual state it may be useful to restore the camera's factory defaults. This returns the capture parameters, image processing, video modes, and image calibration back to the original settings.

To do this, you will need to perform an 'iLoad'. For details on performing an 'iload' follow the 'Step-by-Step Procedures / Live Panel Procedures / Advanced Settings / Performing an iLoad (on a Ph16 Camera)' topic within the supplied PCC Help file.

Back Focus Adjustment

Due to differences in various lenses and other mechanical tolerances it may be necessary to adjust the lens mount on the camera to obtain proper focus. In most cases, changing a lens or replacing a mount will not require adjusting the back focus, though it should be verified. When the focus distance marks on a lens do not match the measured distance between the sensor & subject, it indicates the need to adjust the back focus.

By adjusting the back focus, you are changing the distance of between the flange and the sensor plane. Shims are included to assist in both slight and coarse back focus adjustments.

MAINTENANCE

Sensor Cleaning Technical Tips

- Sensor cleaning should only be attempted by experienced imaging professionals.
- Use a small, but powerful flashlight to look at the sensor and filter, it's easier to spot the dust.
- Sensor cleaning must be done in a clean and controlled environment.
- Always have a forced-air bulb style blower and a clean sensor cloth handy for more serious dirt.
- Sensor cleaning solution (like Eclipse) should only be used if absolutely necessary, and applied with a soft wipe intended for sensor cleaning.
- Cotton swabs aren't a good idea, unless they are wrapped in a sensor wipe material. They leave behind more than they remove.
- Never use your finger to get a speck of dust off, it's going to make things much worse.
- Never blow on the sensor.
- Never put sharp objects near the surface of the sensor.
- Always keep the Phantom body cap on the camera when there is no lens attached.

1 A I N T E N A N C E

CONTACT SUPPORT

GLOBAL HEADQUARTERS

Vision Research, Inc.—Wayne, New Jersey

100 Dey Road Wayne, New Jersey 07470 USA T: +1.973.696.4500

For answers to most questions, please visit us at: *www.phantomhighspeed.com* and search the camera product pages, tutorials, support knowledgebase and FAQs.

SUBMITTING A SUPPORT TICKET

For technical product support, operation and application information or to request an RMA, please submit a ticket by filling out a form at *www.phantomhighspeed-service. force.com* or by emailing us at **phantom-support@ametek.com**.

LIVE CUSTOMER AND TECHNICAL SUPPORT

Serving the Americas:

M-F 8:00 AM to 5:00 PM EST (GMT -4:00) T: +1.973.696.4500 Customer Support, ext. 4002 Technical Support, ext. 4003

Serving Europe, the Middle East and Africa:

M-F 9:00 AM to 6:00 PM GMT +3:00

Vision Research, Inc.—Bucharest, Romania

T: +40 21 210 8587

Serving Asia Pacific:

M-F 8:00 AM to 5:00 PM GMT +8:00

Vision Research, Inc.—Shanghai, China

T: 86-21-58685111, ext. 141

CONTACT SUPPORT

REGULATORY



CE - 2019CE EmissionsCE Compliant - EN 61326-1, Class ACE ImmunityCE Compliant - EN 61326-1, Class AFCCCFR 47, Part 15, Subpart B & ICES-003 Class ASafetyIEC 60950-1



KC - 2022 Country of Origin: USA **R-R-VRi-MiroC321-C321J**: All models Phantom C321 Series KC Emissions KC Compliant - KS C 9832 KC Immunity KC Compliant - KS C 9835

Class A device (Broadcasting Communication Device for Office Use): This device obtained EMC registration for office use (Class A), and may be used in places other than home. Sellers and/or users need to take note of this.

PHANTOM[®] WWW.PHANTOMHIGHSPEED.COM

The contents of this manual are subject to change without notification.

PN: ZDOC-64106-MA-0003 Rev 2 Last Updated: August 2022



100 Dey Road Wayne, NJ 07470 USA +1.973.696.4500